

Product name	X-Safe ImteX	X-Soft ImteX	NeoClean ImteX
Order number			
Materiel	Nitrile	Nitrile	Neoprene
Price per piece	kr.	kr.	kr.
Breakthrough time			
Acetonitrile	1 min	1 min	?
Acetone	?	?	?
Acrylamide, 40%	480 min	480 min	?
Chloroform	0 min	0 min	?
Ethanol, 70%	45 min	?	21 min
Ethanol, 99%	17 min	17 min	?
Ethidumbromide, 5%	480 min	480 min	?
Formaldehyde, 37%	480 min	480 min	121 min
Hydrogen peroxide, 30 %	83 min	280 min	?
Methanol	6 min	5 min	?
Sodium hydroxide, 50 %	460 min	480 min	> 480 min
Sulfuric acid, 98%	7 min	7 min	?

White boxes = not tested for this chemical

**PERMEATION BREAKTHROUGH TIMES ACCORDING TO EN374-3:2003 (MINUTES)**

0	1	2	3	4	5	6
< 10	10-30	30-60	60-120	120-240	240-480	> 480
Not recommended	Splash protection		Medium protection		High protection	

DATA GIVEN IN THE TABLE ABOVE ARE BASED ON RESULTS OF LABORATORY TESTS PERFORMED ON THE PALM AREA OF THE GLOVE OR ARE BASED ON EXTRAPOLATIONS FROM THE RESULTS OF LABORATORY TESTS. THESE TESTS WERE RUN USING STANDARD TEST METHODS THAT MAY NOT ADEQUATELY REPLICATE ANY SPECIFIC CONDITIONS OF END USE. BECAUSE ANSELL HAS NO DETAILED KNOWLEDGE OR CONTROL OVER THE CONDITIONS OF END USE, ANY OF THESE DATA MUST BE ADVISORY ONLY, AND ANSELL MUST DECLINE ANY LIABILITY.